

Installation of Drivers for U90-10 USB Interface

1. Copy the U90-10 folder onto your desktop. You can remove it after installation is complete.
2. Plug in your USB device (U90-10). Windows will open the installation wizard. Using the options shown below, tell windows NOT to search for the drivers. Instead, direct windows to the CDM folder you just placed on your desktop.
3. The installer will run twice, since there are two USB "Ports" one each device.
4. Once installation is complete, the new device will be listed in the System Device Manager as "USB Serial Converter A" and "USB Serial Converter B".

To update drivers for an already-installed device, right-click on each of the two listings in the Device Manager and select "Update Driver..." or "Update Driver Software..." and follow the same procedure as shown below.

WINDOWS 7



WINDOWS XP





